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RUB

GENETIC
PSYCHOLOGY


HOW TO LABEL SAMPLES

a pictured guide

How to label

- use consecutive numbers for identifying your samples, no matter what your subject or patient code is
 - label them on cap and side
- the lab will stick to the numbers you provided when assaying the samples
- make sure to use a waterproof pen that can overcome thawing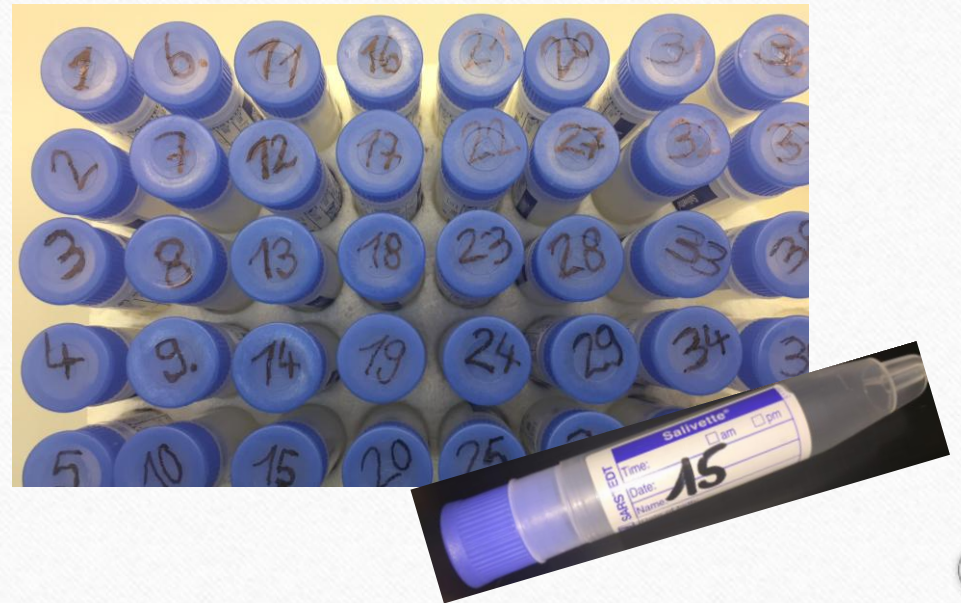

How to pack


- put exactly **20 salivettes** in each zipper bag
- label the bags with a waterproof pen or adhesive label as follows:

bag 1 samples 1 to 20

bag 2 samples 21 to 40


bag 3 samples 41 to 60

to be continued...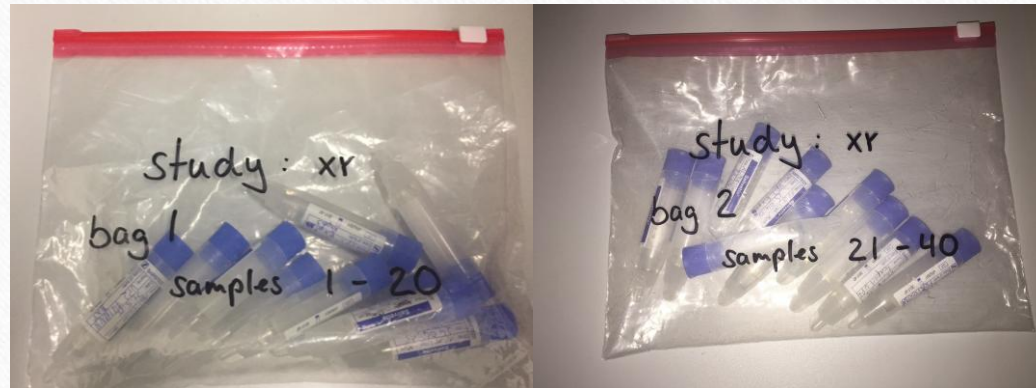

Please do not:

- wrap tape around the tube since it will not fit in the centrifuge anymore
- put the cotton swab in the lower part of the Salivette
 - the swab must be returned to the small beaker/insert
- use green-capped salivettes!
(contains citric acid, interferes with most assays)

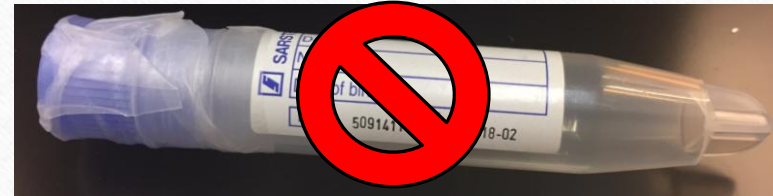How to ship

- Regarding shipping instructions please refer to **"Instructions for the Storage and Shipping of Saliva Samples"**
- please ship the samples to:

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when you need to provide a contact phone number, please use: +49 234 32 24616

Please attach the following statement to your shipment:

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
Q&A

How soon will I
receive my results?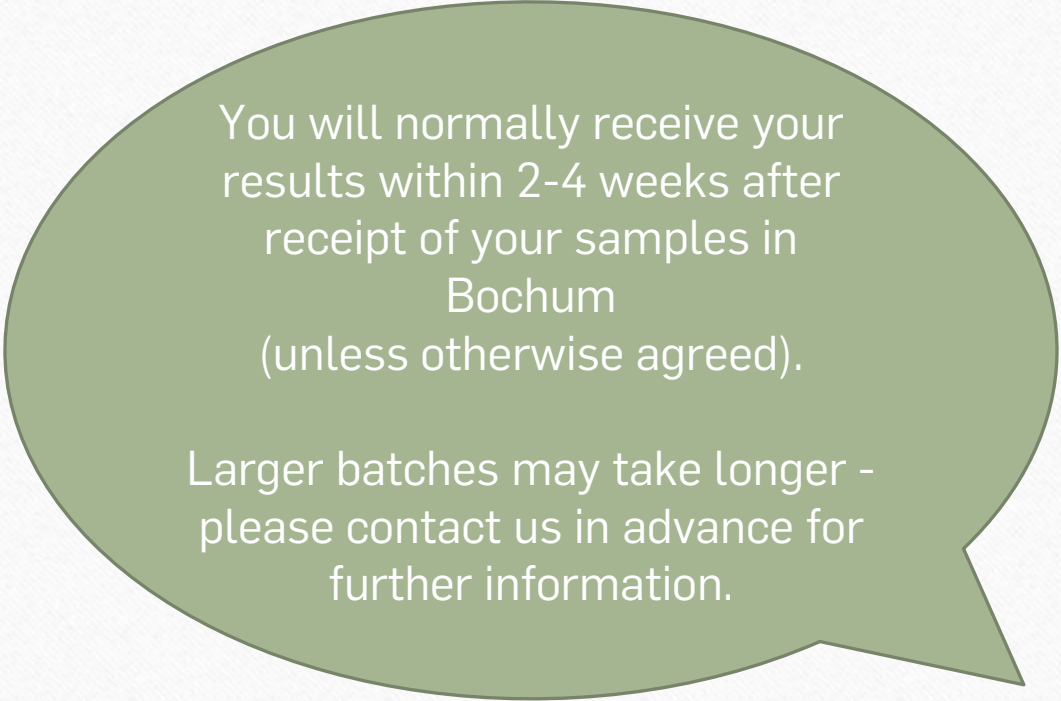


You will normally receive your
results within 2-4 weeks after
receipt of your samples in
Bochum
(unless otherwise agreed).

Larger batches may take longer -
please contact us in advance for
further information.

Q&A

How do I ship my samples to you?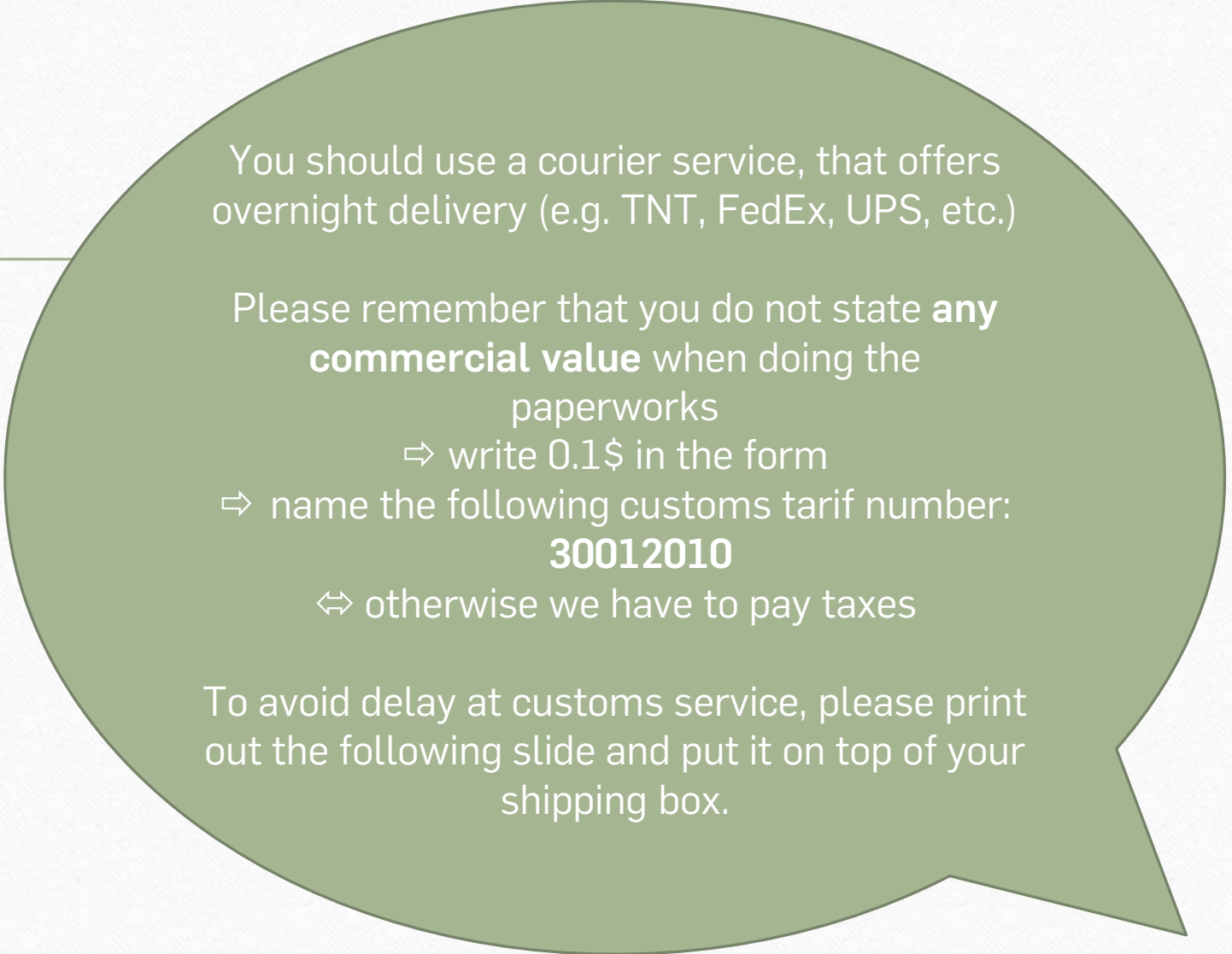

You should use a courier service, that offers overnight delivery (e.g. TNT, FedEx, UPS, etc.)

Please remember that you do not state **any commercial value** when doing the paperworks

⇒ write 0.1\$ in the form

⇒ name the following customs tariff number:

30012010

⇔ otherwise we have to pay taxes

To avoid delay at customs service, please print out the following slide and put it on top of your shipping box.